ELBER

M E D I C A L

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ELBER

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26	MAINTENANCE AND PREVENTIVE VERIFICATION.....	39
27	EMERGENCY SYSTEM (see availability of optional)	39
28	PROBLEMS AND PROBABLE CAUSES.....	41
29	PRODUCT DISPOSAL	41
30	WARRANTY	41
31	PROCEDURE OF HANDLING THE DIAL TELE ALARM 6 NUMBERS (<i>see availability of the optional</i>).....	42
31.1	Register phone numbers	42
31.2	Erase numbers for memory	43
31.3	Test tele alarm memories.....	43
32	PROCEDURE OF HANDLING THE DIAL TELE ALARM 9 NUMBERS (<i>see availability of the optional</i>).....	44
32.1	Programming a phone number	44
32.2	Programming a phone number with pause 2 sec. (for calls via PABX)	44
32.3	Programming a phone number for call via economy line.....	44
32.4	Programming a phone number for long distance calling.....	45
32.5	Deleting a phone number	45
32.6	Testing phone numbers.....	45
32.7	Changing the installer password	45
32.8	Dialing process	45
32.9	Dial-in interrupt modes	45

32.10	TOTAL RESET (All dialer parameters return to Factory Standard).....	46
32.11	Audible signaling (Buzzer).....	46
33	GSM DIALER HANDLING PROCEDURE (see availability of optional).	46
33.1	Installing the SIM CARD	47
33.2	Signs	47
33.3	Programming the GSM Dialer	47
34	WIFI MODULE HANDLING PROCEDURE (<i>see availability of optional</i>).	48
34.1	Set up the Wi-Fi module	48
34.2	Historical.....	49
34.3	Settings.....	50
34.4	Register mobile phones and emails.....	50
35	SURGE PROTECTION (DPS).	50
35.1	DPS operation	51
35.2	Connection and attachment of the DPS ELBER kit	51
36	INVERSÃO DA ABERTURA DA PORTA (ver disponibilidade do opcional).....	53
37	CARACTERÍSTICAS DE CONSERVADOR.....	55
37.1	Embalaje	55
37.2	Transporte	56
37.3	Almacenamiento	56
37.4	Manipulación	56

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DEAR USER!

You have just acquired a conservative for the storage of immunobiologicals, blood products, medicines, thermolabes and the like. This product was developed with ELBER technology and functionality. We are grateful for your choice.

It is essential to carefully read the instructions contained in this manual.

Following these guidelines, you will be able to know better the operation of the product, enjoying its benefits and prolonging its durability. The instruction manual must always be accessible near the product, so that any questions can be answered immediately.

ANVISA Registration: 80698750001 and 80698750002.

Authorization (AFE) ANVISA: 8069875

19 CHARACTERISTICS OF CONSERVATIVE

Elber conservatives were developed to store immunobiologicals, blood products, drugs, thermolabes and similar to a stable, homogeneous and uninterrupted temperature. It is indicated for hospitals, pharmacies, health clinics, vaccine clinics, blood banks and laboratories in general.

Works through electric power 110-127 VAC or 220-240 VAC.

ELBER conservatives are designed to ensure that the internal temperature of the chamber is always stable and homogeneous. Equipped with a security system, the conservative Elber remain functioning even with a lack of commercial electricity (see availability of the optional). Elber conservatives have a microprocessor electronic controller that controls the internal temperature of the camera (cooling as needed) and informs the user of all events through audible and visual alarms and electronic records.

Elber conservatives can be equipped with optional amusements, see in the technical specification sheet, which accompanies your conservative, all product characteristics.

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After unpacking the product, any material in the package should be sent for recycling or further reuse.

19.2 Transport

Elber sums must be transported upright, maintaining packaging conditions according to the label.

19.3 Storage

ELBER conservatives should be stored in clean, dry, sun-protected environments and excessive humidity.

The conservative should not be stored for more than two months without being used. If not used, the battery will discharge.

19.4 Manipulation

ELBER conservatives should only be handled by trained professionals who fully meet the specifications of the product instruction manual.

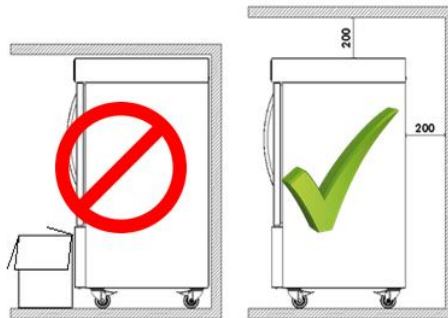
20 INSTALLATION AND FIXATION OF THE CONSERVATIVE

The conservative should be positioned on a flat floor and leveled. The feet have casters that allow you to position it easily. To move the conservative the casters must be unlocked and locked when the conservative is positioned. Moving the conservative with the locked casters may damage the casters and the floor.

For large conservatives (1000, 1200 and 1500 liters), after positioning the conservative, level the product using the leveling feet of the sides (near the casters). The central supports should be threaded until you touch the floor, do not use these as levelers, only as supports.

It should be installed preferably protected from the sun and in a place where there is air circulation. Very small rooms should have openings for air exchange.

We do not recommend installing the conservative near heat sources (heaters and greenhouses). In the ventilation grilles of the conservative should there must be space of at least 20 centimeters from the wall or any obstacle, to allow the ventilation of the compressor unit and the electronic system. For conservatives equipped with emergency system of 72h, the machine base is larger and has ventilation at the top and rear side, in this case the conservative can be leaned against the wall at the back.

21 ELETRICAL INSTALLATION

The conservative is prepared to run connected on 110-127 VAC or 220-240 VAC power.

Before turning on the battery, check that the voltage indicated on the product coincides with the voltage of the mains

The power cord must be plugged into a permanent fixed outlet on the wall or worktop in accordance with current regulations. Do not use extensions or adapters..

Note the steps below to turn on the conservative correctly.

1º - The voltage 110-127 VAC or 220-240 VAC is factory adjusted;

2° - Plug the plug into the socket - the conservative will not work yet, as the general switch is missing;

3° - Turn on the general switch of the conservative;

4° - When turning on the conservative, the **Sound Alarm** will sound until the temperature inside the camera is within the ranges set on the digital controller.

You can cancel the audible alarm for a predetermined time by pressing the mute key (ELV5 and touch screen) or any key (ELV4) of the controller. When the alarm is sounding and is turned off, it will sound again after the set time if the temperature remains outside the programmed range. The alarm will sound until the conservative reaches the normal operating range (on average 2 hours without charge).

To turn off the conservative, when equipped with emergency system, it is necessary to turn off the general switch. It is not enough to remove the plug from the socket, as the conservative will continue to work by the battery (*see availability of the optional*).

The use of the Surge Protection Device (DPS) SPW 275-45 is indispensable. This device protects the conservative from surges and tensions, preventing the burning of it.

Attention

This conservative should be connected in a ground to prevent damage to the product and the user. Ground wire connection is required at any voltage (110-124 VAC or 220-240 VAC). The lack of grounding and the use of the DPS device compromises the coverage of the conservative warranty. To perform the grounding and electrical installation, as well as the DPS, follow the standards of ABNT-NBR 5410, or call a responsible professional.

22 OPERATION

This conservative is prepared for the conservation of immunobiologicals, blood products, medicines, thermolabes and the like.

It guarantees a constant temperature at any time of the year and works even when there is a lack of electricity (see availability of the optional).

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The **guarantee of a uniform temperature** is made by the digital controller that maintains what has been programmed. The controller is programmed to keep the camera at a temperature between +2.0°C and +8.0°C, both on cold and hot days. At any time of the year, winter or summer, the conservatory will automatically adjust.

The temperature will always remain in a defined range, ensuring the best properties of the stored material.

23 AUDIBLE ALARM

For convenience, the conservative has an audible alarm that is triggered when the temperature is outside the programmed limits or other anomalies. This can occur if the door is open too long or by some malfunction. So that in a timely manner some action can be taken so that the products in storage are not lost. The alarm will be activated only after the predetermined time with temperature out of schedule.

Function delay alarm: When turning on the conservative alarm will sound. To suspend the alarm for the predetermined time, press the mute key (ELV5 and touch screen) or any key (ELV4) of the digital controller. Repeat the operation as many times as necessary until the conservative reaches the programmed track, where it will no longer fire. To change the time between shots, see regular time function to trigger the alarm.

24 BEST USE OF CONSERVATIVE AND RECOMMENDATIONS

- Avoid leaving the door open. D
- In use, always keep products inside the conservative, even if they are bottles with water. This load of products helps to keep the temperature stable inside the conservative. D
- Before placing the vaccine in the conservative check that the operation is correct and the air circulation fan on the inner bottom is working. D

- Avoid leaving the conservative door open for too long or open it constantly;
- Keep the conservative running for at least 48 hours before placing products in the conservative, to check that the general operation is correct, that the internal fan is circulating the air and that the temperature is within the set parameters;D
- In use, always keep products inside the conservative, even if they are bottles with water. This load of products inside helps to keep stable the temperature inside the conservative;
- Internal ventilation flaps must not be obstructed, constantly check whether stored products are obstructing the fins and, consequently, internal ventilation of the product;
- Shelves or drawers must not be lined with plastics, towels or the like, which compromises the internal ventilation of the product;
- On days with a lot of moisture, there may be condensation of air moisture (sweat) on the door or at the outer edges of the conservative (near the gasket). In places with room temperature controlled by air conditioning, this feature is practically extinct.
- For conservatives equipped with emergency system, do not perform battery life tests without technical monitoring Elber, face-to-face or online. Only technicians can guarantee the recording of the results, disable the messaging system and ensure that the battery does not suffer damage (see availability of the optional).
- Only trained persons should operate the product.

25 DEFROSTING AND CLEANING

Elber snowpacks are equipped with dry automatic defrosting. To clean the conservative, first turn off the GENERAL SWITCH. It is not enough to unplug it, as it will continue to operate by battery (see availability of the optional).

Use damp cloths with water and neutral detergent. Do not use splashes or water hoses, excess water may compromise the operation of the product.

Do not use chemicals (alcohol, solvents, etc.) or abrasive products and hard sponges. Non-recommended procedures may stain and damage the surfaces and coating of the refrigerator.

During cleaning, remove all materials stored in the conservative and place them in other equipment or thermal box with cryogenic material.

Important! After a period of conservative inactivity (for cleaning, for example), when turning on, the conservative can reach temperatures below the beloved program the electronic controller. Therefore, whenever you turn off the conservative, after reconnecting, wait at least 2 hours to put load on the conservative again. After 2 hours, clear the controller data and start normal use.

26 MAINTENANCE AND PREVENTIVE VERIFICATION

- From 12 months on it is important to check for soot or dirt that may be clogging the condenser;
- After 03 years of use of the equipment, it is recommended to test the battery life of the emergency system (test consists of removing the equipment from the outlet for 2 hours, it must be monitored, to see if there is oscillation of internal temperature of the equipment) Periodicity - every 6 months;
- After 05 years of use of the equipment, it is recommended to perform the replacement of the internal fans and fan of the condenser of the chamber;
- According to the special characteristics of the equipment and/or customer's requirements, it can be adapted preventive maintenance system according to customer's requirements;

27 EMERGENCY SYSTEM (see availability of optional)

When equipped with emergency system, in lack of conventional electric power 110-127V or 220-240V, the elber conservatives continue to operate (see autonomy in the sheet of technical specifications of the product), keeping all its electronic and cooling functions in operation.

The activation of the emergency system is automatic. The autonomy of the emergency system varies for each model, can be increased or decreased, as needed by the customer.

Elber's conservatives have maintenance-free sealed batteries with an automatic charger.

Battery recharge time varies between 10 and 20 hours.

Important: Battery life depends on the temperature conditions of the conservative installation site, the number of times the equipment door is opened with the emergency system in action (without commercial power), and several other factors.

The emergency system is an optional, check availability of the optional.

Attention: In new conservatives with emergency system, in the first hours after the restoration of commercial electricity, the conservative can exude a chemical odor, caused by a slight heating in the battery charger. This condition is normal and only occurs in the first cycles of operation of the emergency system (after the lack of commercial electricity). In this condition, it is recommended that the environment where the conservative is installed be opened until the odor is eliminated.

28 PROBLEMS AND PROBABLE CAUSES

FAULT	SOLUTION
Conservative doesn't care	<ul style="list-style-type: none"> - Check that the voltage of the commercial electricity network is the same as the conservative one; - Check for commercial electricity; - Check that the general key is on; - Check that the plug is plugged into the outlet; - Check that the protective fuse is blown; - Check for any warnings on the digital controller.
Conservative does not cool;	<ul style="list-style-type: none"> - Check that the compressor is on (place your hand on the compressor and check for vibration and light heating); - Check that the internal ventilation is working; - Check the door closing (door gasket cannot have gaps in relation to the cabinet); - Check for any warnings on the digital controller.
Conservative alarming by high temperature	<ul style="list-style-type: none"> - Check that a large product load has been loaded less than 2 hours (the conservative can take until it reaches the ideal temperature when a large amount of product is placed at once); - Check the door closing (door gasket must not have gaps in relation to the cabinet). - Check for any warnings on the digital controller.
Internal temperature of the varying conservative.	<ul style="list-style-type: none"> - Check that the internal fans are working; - Check that the internal ventilation fins are free (they must not be obstructed);

29 PRODUCT DISPOSAL

Out-of-use or abandoned refrigerators can be dangerous, especially for children, as they can get trapped inside, at risk of shortness of breath. Before disposing of your old refrigerator cut the power cord and remove the door(s). Leave the shelves in place so the children can't get in.

In case of battery disposal of the conservative, forward the discarded battery to a nearest pick-up point.

In case of disposal of the product, all materials must be sent for recycling.

30 WARRANTY

Conservative warranty is 1 year.

Attention: The battery warranty term is next to it, and cannot be lost as it contains the serial number that matches the battery number (see availability of the optional).

31 PROCEDURE OF HANDLING THE DIAL TELE ALARM 6 NUMBERS *(see availability of the optional).*

The tele alarm works through a telephone line. Makes calls to the numbers recorded when the temperature drops out, or if the battery voltage is low. For the dialer to work, you need to connect a telephone cable.

31.1 Register phone numbers

The tele alarm has memory to record up to 6 phone numbers, with up to 16 digits each.

To register a phone number, press the "PROG" key to enter programming mode. Type "1", which corresponds to the desired function, and then press the memory number the phone number should occupy, from 1 to 6. Enter the phone number and press "ENTER" to confirm. Press "ENTER" again to exit programming mode. For example:

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When the line has a PABX extension, press "0" to pick up the outside line, and the "PROG" key, to insert a 2-second pause before the phone number. Example:

If a number is written to an already occupied memory, the new number replaces the previous one, which will be deleted automatically.

We recommend recording numbers from at least two different carriers, if the carrier fails and it is not possible to perform operation for one, the dialer can complete the connection to another.

In case of firing the tele alarm, it will call all numbers, according to the registration sequence. The dialer will continue to sequence calls as long as there are shots.

31.2 Erase numbers for memory

Press the "PROG" key to enter programming mode. Type "2", which corresponds to the desired function, and then press the memory number that corresponds to the phone number you want to delete (from 1 to 6).

Press "ENTER" to confirm, and again to exit programming mode. Example:

31.3 Test tele alarm memories

Press the "ENTER" key and then the memory number corresponding to the phone number you want to test. The LED will flash twice, rise and start the test call. If the memory is empty, the LED rises by 1 second and the test will be finalized. Ex.:

To cancel the test, press the "PROG" key.

THE OTHER TELE ALARM FUNCTIONS ARE ALREADY FACTORY PROGRAMMED, NOT BEING NECESSARY TO CHANGE.

32 PROCEDURE OF HANDLING THE DIAL TELE ALARM 9 NUMBERS *(see availability of the optional).*

The tele alarm works through a telephone line. Makes calls to the numbers recorded when the temperature drops out, or if the battery voltage is low. For the dialer to work, you need to connect a telephone cable.

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32.2 Programming a phone number with pause 2 sec. (for calls via PABX)

Press # + Installer Password + Index (01 to 09) + 0 + E + Phone Number + # + #.

If you need to use the characters * (asterisk) or # (cerquilha), proceeded in the same way, pressing, respectively, the P or F key in the position where they should be inserted.

32.3 Programming a phone number for call via economy line

For calls of this type, where it is necessary to type *15 seconds per pause of 6 seconds before completing the dial, press # + Installer Password + Index (01 to 09) + P + 15 + E + E + E + Phone Number + # + #.

Phone numbers older than 22 digits are not allowed. When the 22nd digit is entered, the system automatically writes its value, and it does not have to press # to store it.

32.4 Programming a phone number for long distance calling

Press # + Installer Password + Index (01 to 09) + 0 + Carrier Number + Phone Number + # + #.

32.5 Deleting a phone number

Press # + Installer Password + Number Position you want to delete (Index 01 to 09) + * + # + # #

Another way to delete a programmed number is to access the desired index and omit the value of the phone number by pressing #. Be careful when accessing phone indexes (01 to 09) accidentally, as you will need to re-enter their value.

32.6 Testing phone numbers

With the phone line properly connected and the dialer in user mode (off-schedule), press F + Desired Memory Position (1 to 9 - without digit 0).

The dialer will make a call to the stored number, playing a siren sound, on the phone for 60 seconds. During this time, the LINE LED will remain lit. To stop a test in progress, press any key.

32.7 Changing the installer password

The installer password is used to access The Programming Mode and its factory value is 4321. To change, press # + Installer Password (current) + 10 (index) + New Installer Password (4 digits) + #.

32.8 Dialing process

The dialer runs through a call cycle that starts with the phone registered in the first memory position (index 01) and ends with the phone registered in the last memory position (index 09), performing up to 2 dial cycles.

Ex: Start of the first cycle / MEM1... MEM2... MEM3... MEM4... MEM5... MEM6... MEM7... MEM8... MEM9... / End of first cycle / MEM1... MEM2... MEM3... MEM4... MEM5... MEM6... MEM7... MEM8... MEM9... / End of the second cycle and the dialing process.

32.9 Dial-in interrupt modes

Simultaneous Mode: The dialer will make calls only as long as the alarm or fence is set off.

Persistent Mode: The dialer will make calls even if the trigger ceases to occur, completing the 2 cycles.

Press # + Installer Password + 12 (index) + 1 (Simultaneous Mode) or 2 (Persistent Mode) + #.

32.10 TOTAL RESET (All dialer parameters return to Factory Standard)

Turn off the dialer power and hold down keys 7 and 9. Realistically, temporarily dial the dialer until you hear 3 short beeps. Release keys 7 and 9 and realistically dial.

32.11 Audible signaling (Buzzer)

1 Long Beep = Error Tone (incorrect command or procedure).

1 Short beep = Indicates the pressing of a key.

3 Short beeps = Confirmation Tone (command or procedure accepted).

THE OTHER TELE ALARM FUNCTIONS ARE ALREADY FACTORY PROGRAMMED, NOT BEING NECESSARY TO CHANGE.

33 GSM DIALER HANDLING PROCEDURE (see availability of optional).

The conservative Elber can be equipped with GSM dialer, which sends SMS messages and makes calls in case of events. Below are the main instructions related to this item.

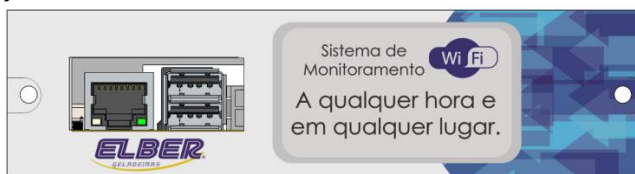
- In case of long-distance connections, it is mandatory to choose the operator and the area code, or else the call will not be completed.
- If a landline is registered, it will only receive trigger alert and or low battery.

To delete registered phone numbers, enter the "02" function, select the memory you want to erase (1 to 6) and press "E" to confirm.

To test the memorized numbers, press the "E" key and then the memory number (1 to 6). If there is a number stored, the "E" key will remain lit, then start sending the SMS with bitonal alert link. If the memory is empty, an error BIP is issued and the test is terminated.

For more detailed information on GSM dialer operation, refer to the manufacturer's manual www.sulton.com.br.

34 WIFI MODULE HANDLING PROCEDURE *(see availability of optional).*

The Wi-Fi module allows the monitoring of the system via the Internet. Allows you to obtain reports, system information and register mobile numbers and emails to be notified in case of problems, all through a computer with internet access.

34.1 Set up the Wi-Fi module

To configure the module, you will need a pen-drive and a computer. Follow the steps below:

1. Insert the pen-drive into your computer and create a notepad file with the name *internet.txt*. This file cannot be saved within a pen-drive folder.
2. Type in the first line the name of the Wi-Fi network that the module must be connected to, followed by a comma and password. Example: *redeelber, senhaelber*.

Note: Do not enter space after the comma.

3. Remove the pen-drive from the computer and insert it into one of the Wi-Fi modules ports.
4. Wait for the ELV5 controller display to inform you that the network has been set up and that the drive pen-drive can be removed.
5. The module will start sending information for remote monitoring. To gain access to online data, send an email to at04@elber.ind.br and at05@elber.ind.br requesting registration.

After accessing the account, the screen below will appear, in addition to the temperature, voltage and GSM signal level graphs, when available.

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To register numbers and emails and generate the reports, follow the instructions below.

Attention!

When possible, use network cable on the WiFi system.

NOTE: Monitoring system that depends on transmission and or storage of data by Wi-Fi, GPS, SMS, WhatsApp, or satellite are subject to change of technologies, licensing and possible related costs, for maintenance of monitoring.

34.2 Historical

By clicking *on history*, you can get the reports of the temperature, battery voltage and electrical power source charts. After clicking, select the desired period and click *filter*. Charts and data table will appear at the end of the page. If you want, click Print.

34.3 Settings

The settings should be changed only by the company Elber.

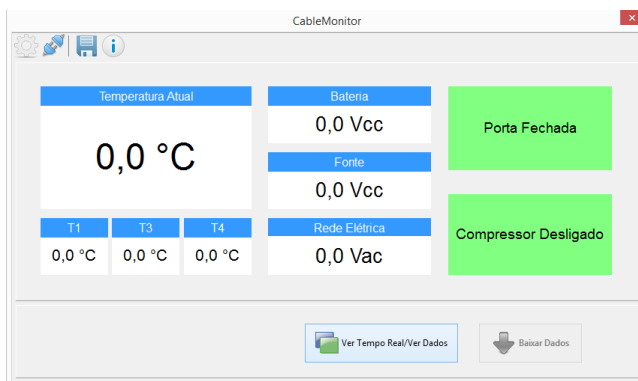
34.4 Register mobile phones and emails

By clicking *on email*, you can register emails to receive messages informing problems, such as temperature incurring and lack of electricity. After clicking email, click add and enter as many emails as you want.

By clicking *on mobile phones*, you can register mobile numbers to receive messages informing you of problems. After clicking mobile phones, click add and enter as many numbers as you want.

INTERFACE MODULE *(see availability of the optional)*.

The purpose of this function is to download the data straight to the computer without the need for a thumb drive and monitor conservative events in real time. With the computer plugged into the conservative via USB cable, you must use the software "Cable Monitor". After clicking on "connect", the software will show the real-time events of the conservative. To download the report, the user must click "Download Data". When installing The ELV5 CHART is automatically downloaded and installed the software "Cable Monitor".

35 SURGE PROTECTION (DPS).

DPS Elber is an outbreak protection device (DPS). It is an equipment that aims to detect the surges in the electrical network and divert the surge currents, it

is an indispensable device for the use of Elber products. It can be installed in the circuit breaker box on site or via the DPS ELBER Kit.

Maximum discharge current of the device: 45 kA

35.1 DPS operation

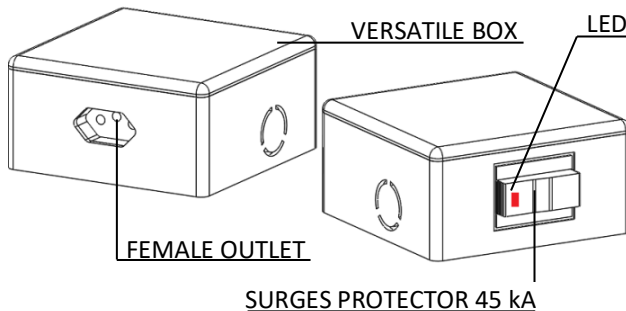
The DPS Elber is connected to the electrical network and responsible for the protection of outbreaks, when subjected to an "Electrical Surge" the DPS will stabilize the electrical current that the equipment will receive.

WARNING: The Surge Protector when subjected to a very large power variation may be damaged, it will signal through an LED that will light up, and replacement of the protector will be required.

WARNING: To perform the DPS installation follow the standards of ABNT-NBR 5410:2004, DPS section, or activate a responsible and qualified professional.

35.2 Connection and attachment of the DPS ELBER kit

For connection of the DPS Kit, it is essential to shut down the electric current!


CONNECTION OF THE DPS ELBER KIT:

The DPS Elber kit consists of an Surge Protector, a female socket for connecting the equipment to be connected and a box.

The DPS Elber kit must be connected directly to the mains, respecting all required safety standards. It is indispensable to connect the "Earth" to the kit, so it can provide all the safety proposed to your equipment.

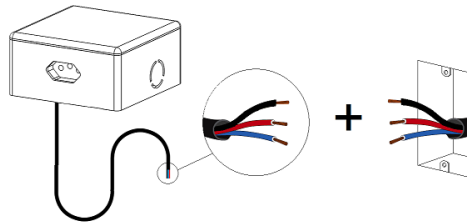
The kit comes with a Phase, Neutral and Earth connection cable, which will exit from the bottom of the DPS kit.

WARNING: The correct insulation of the wires is indispensable, either by an insulating tape or retractable thermo spaghetti.

DPS KIT FIXING:

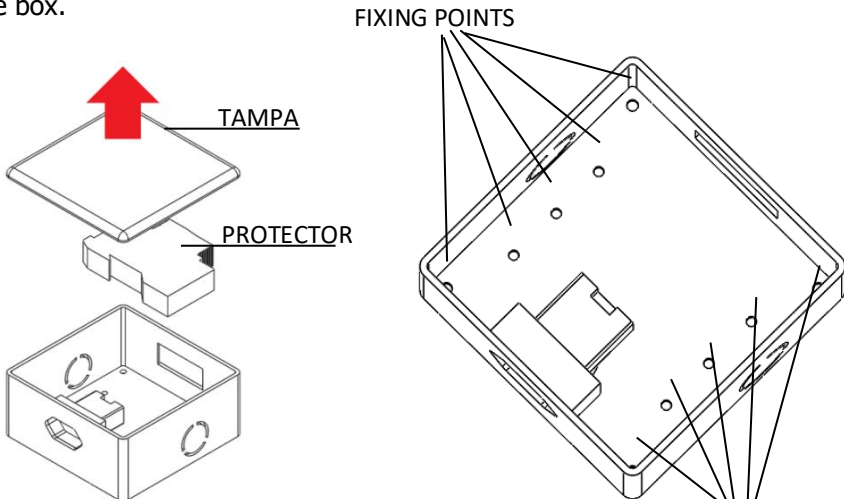After connecting the DPS to the mains, and still with the electric current off, remove the top cover of the Kit box, if necessary, also remove the Surge Protector, for better access to the attachment points located at the bottom of the box.


To secure the DPS Kit housing, simply position it to the desired location and insert the screws through the attachment points.

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The conservative Elber, have the technology and practicality that allows the inversion of the opening of their doors. To do this, follow the steps below:

1. Remove the top cover of the conservative, to gain access to the top hinge of the door;
2. Tilt the front cover of the base to gain access to the inner cover of the base;
3. Remove the inner cover of the base to gain access to the lower hinge;
4. Dispatch the electrical connection of the door with the base of the conservative;
5. Remove the lower and upper hinge (care is required in this process, as the door must be held so that it does not suffer any kind of damage);
6. Rotate the 180° port for its installation so that its opening is reversed;
7. Install the hinges in reverse, and the upper one will become the lower and the lower will become the upper hinge;
8. After fixing the hinges, reinstall the door to the conservative;
9. Re-place the electrical connection of the door to the base of the conservative;
10. Reassemble the base covers, as well as the top cover of the ceiling;

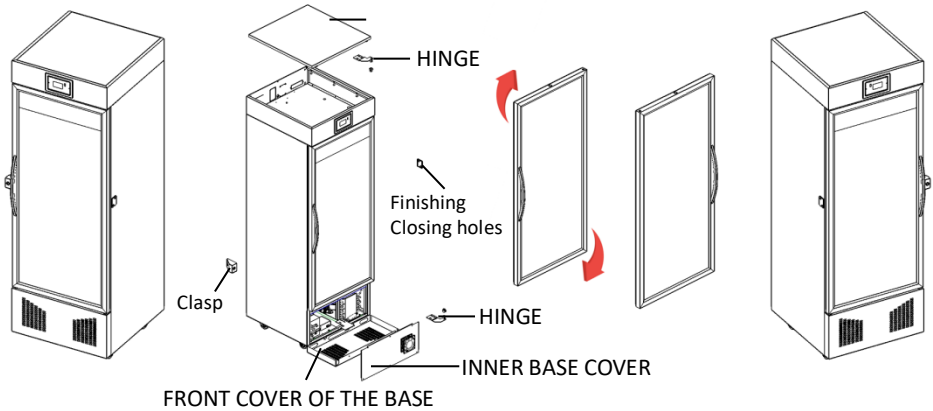CEILING COVER

(*) For conservatives who have a lock on the door, they will be prepared to meet the inversion of the door opening as well, just remove and install the clasp and finish the holes on the other side of the conservative.

Next to the conservative, a new adhesive of the door will be sent in which it can be applied after the inversion of the same and removal of the previous adhesive.

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